

Advanced Cell Culture Solutions for Your Seahorse XF Analyzer

Agilent sensor cartridges and cell culture microplates





Gain a Clearer Picture of What Your Cells Are Doing

Comprising instruments, assay kits, software, and consumables, Agilent Seahorse XF platforms are designed to support you from sample to insight.







Sensor cartridges for reliable kinetic measurements

Cellular oxygen consumption (respiration) and proton excretion (glycolysis) cause rapid changes to the concentrations of dissolved oxygen and free protons. Agilent sensor cartridges for Seahorse XF analyzers ensure sensitive, kinetic measurements to simultaneously analyze both metabolic energy pathways.

Key features

- Patented transient microchamber improves analytical sensitivity and signal-to-noise of OCR and ECAR (PER and gPER) readings.
- Solid-state sensor probes containing polymer-embedded fluorophores detect oxygen and proton levels in cell culture media. The probes reside 200µm* above the cells and take measurements every few seconds.
- Integrated injection ports—a patented design—let you add compounds, stimulators, inhibitors, substrates, and beads conjugated with antigens.
- Sensors are spotted onto the cartridge using both manual and machine vision inspection.
- Every lot is 100% quality controlled for performance and lot-to-lot consistency.

* Microchamber volume differs by tissue culture plate type.

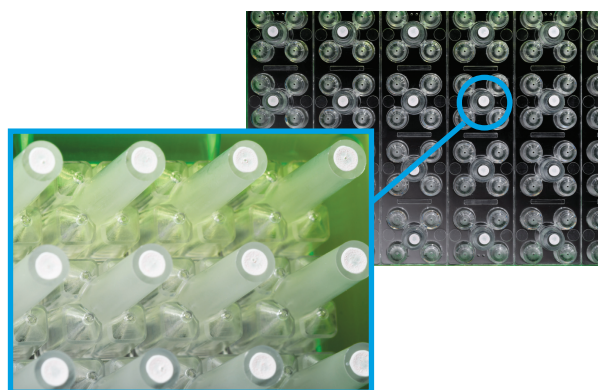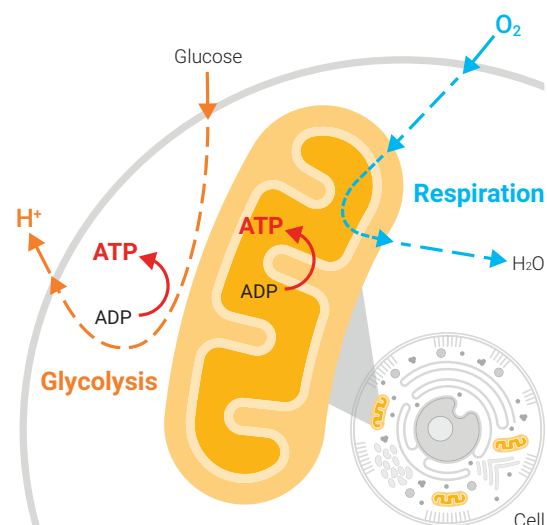


Agilent Seahorse XF FluxPaks Take You from Sample to Insight

Convenient fluxPaks contain the essential consumables you need to perform XF assays.

Sensor cartridges

The essential component for simultaneously measuring bioenergetic pathways.

- Non-invasive sensors to measure dissolved oxygen and free protons
- Four injection ports
- Lot-to-lot consistency
- High-precision manufacturing process
- One-year shelf life from manufacturing date
- Patented design

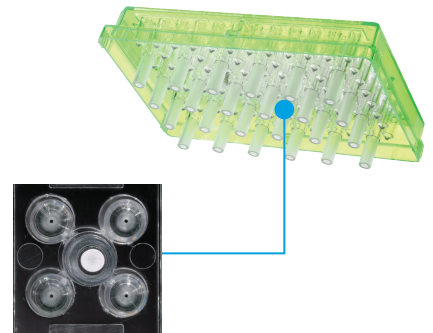

Cell culture microplates


Choose from many plates to suit your application and sample type. See page 8 for a complete list.

- Tissue culture treated and irradiated for sterility
- Tapered walls to facilitate cartridge alignment, especially at low volumes
- Well “bumps” define microchamber height and volume
- Conforms to ANSI/SLAS microplate standards for footprint dimensions and well positions.
- Polystyrene material
- Living cells, isolated mitochondria, islets, and spheroids
- Poly-D-Lysine 96- and 8-well formats



Calibrant and calibration plates

- Premixed calibrant: ready to use and optimized for XF sensor cartridge calibration
- Calibration plates: suitable for calibration and hydration
- Hydrobooster: Optimize hydration workflow

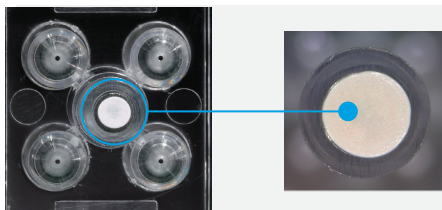


Precise sensor cartridge spotting

High-resolution machine vision imaging inspects the spot quality on every cartridge.

Seahorse XF Technology: A Closer Look

The integrated drug delivery system lets you sequentially add up to four compounds per well, so you can perform a complete mitochondrial or glycolytic kinetic assessment.

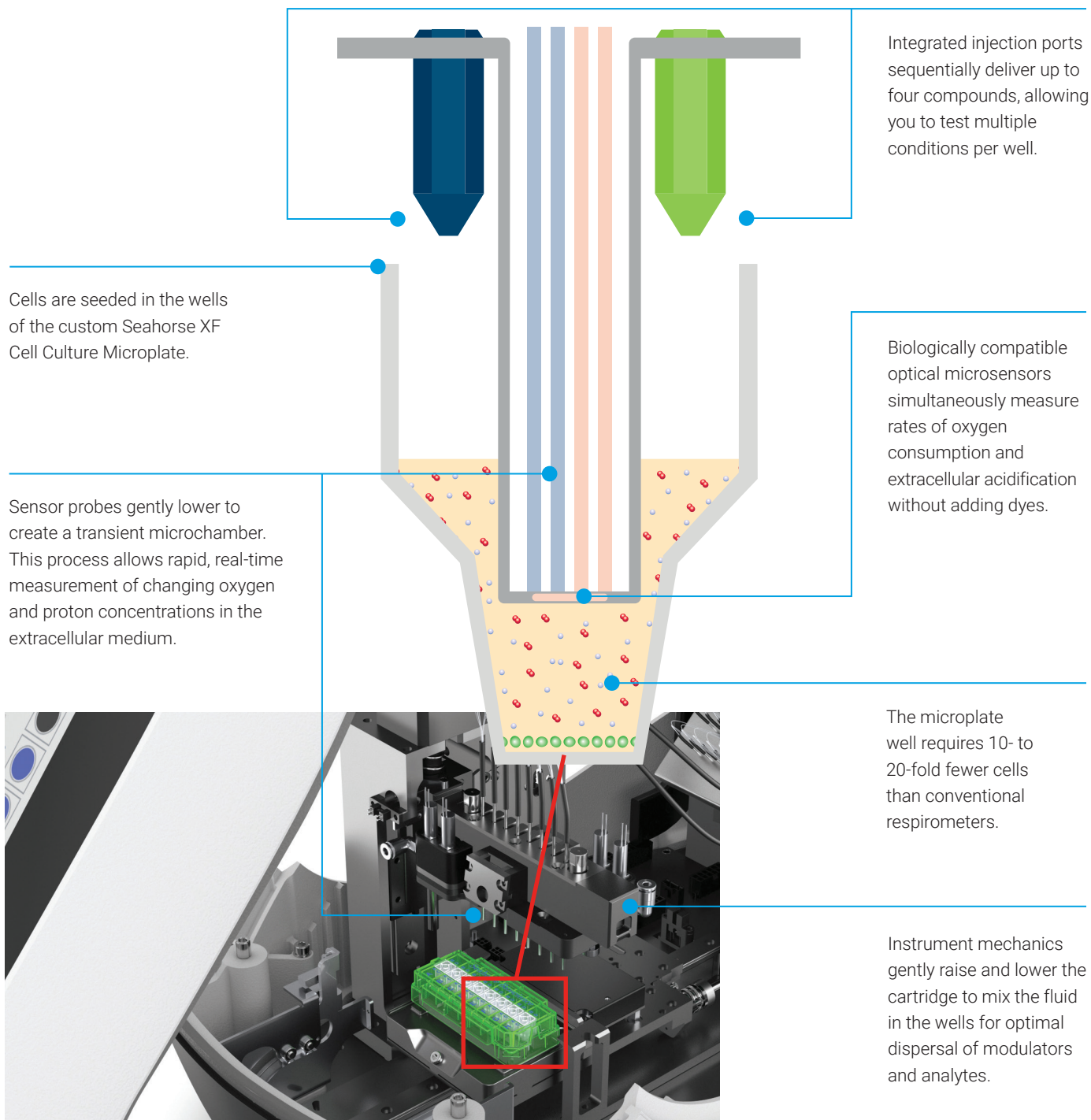


What happens during a measurement cycle?

Each measurement cycle amounts to one time point; over time, the kinetic data is plotted and displayed.


Mix

When the assay starts, the sensor cartridge probes lower into each well and repeatedly mixes the culture medium.

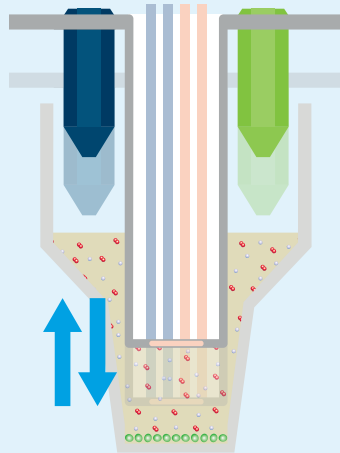


Measure

The probes lower to form the extremely small volume of the transient microchamber.

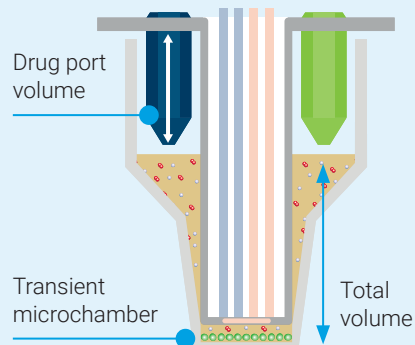

Inject


After a series of basal measurements, constituents are injected from the drug ports into the culture medium at user-programmable intervals.

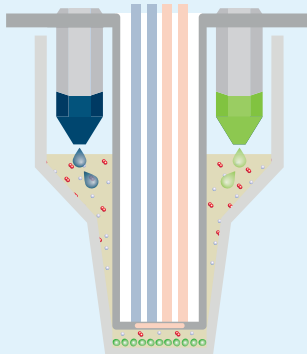


Cell Culture Microplate Types and Specifications




Standard Microplate

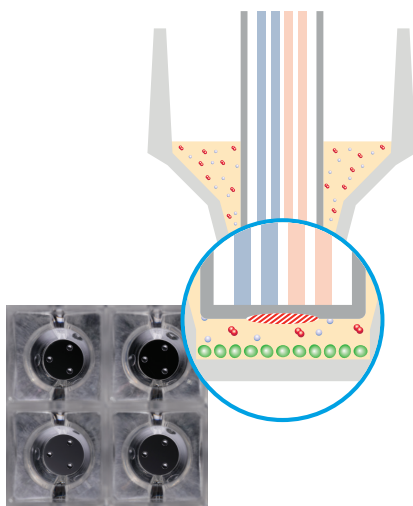


Islet Capture Microplate

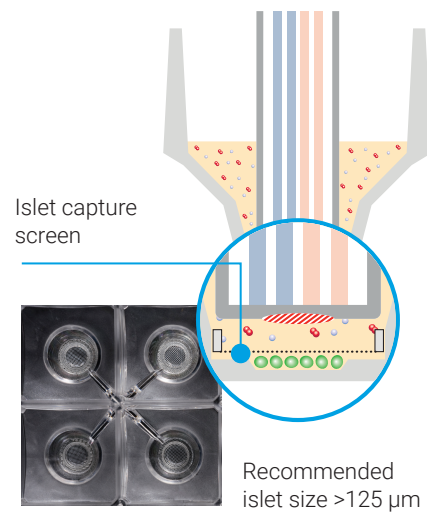


Spheroid Microplate

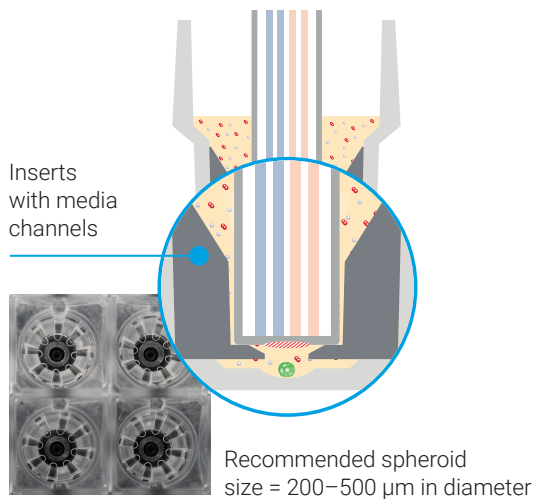

XF HS Miniplate

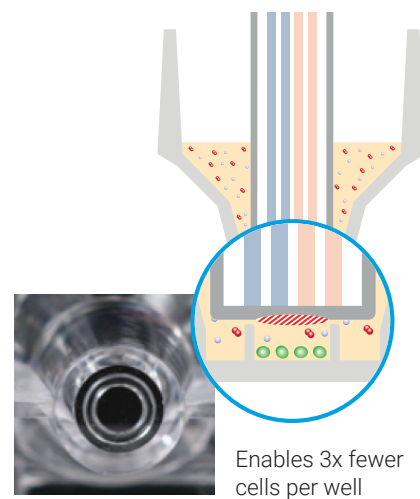

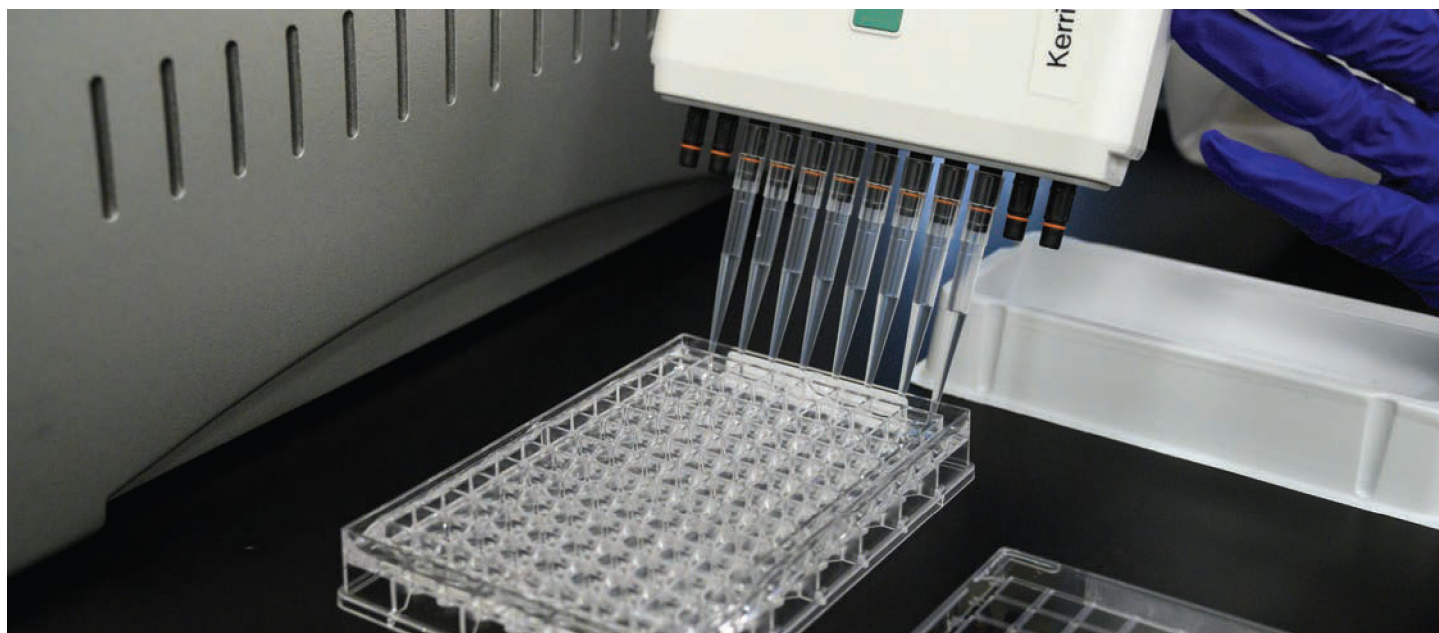

Seahorse XF Cell Culture Microplates and Miniplates can be modified with substrates to improve cellular adherence. You can purchase precoated Poly-D-Lysine plates or modify manually (such as collagen or Cell-Tak).

Analyzer	Plate Type	Number of Wells	Total Volume (μL)	Microchamber Volume (μL)	Height of Microchamber (μm)	Material	Cell Types	Isolated Mitochondria	Part Number
XF96, XF96 or XF Pro	V3 PS	96	250	2.28	200	PS	2-D, monolayer	Yes	101085-004
XF96 or XF Pro	Spheroid	96	250	1.88	850	PS	3-D, spheroids	No	102978-100
XF96, XF96 or XF Pro	V3 PS (+PDL)	96	250	2.28	200	Poly-D-Lysine on PS	2-D, suspension and monolayer	Yes	103729-100
XFe24 or XF24	V7 PS	24	1000	5.65	250	PS	2-D, monolayer	Yes	100777-004
XFe24 or XF24	V28 PS	24	1000	5.65	800	PS	2-D, monolayer	No	100882-004
XFe24 or XF24	Islet	24	1000	16.6	250	PS	3-D, islets	No	101122-100
XF HS Mini or XFp	V3 PS	8	250	2.28	200	PS	2-D, monolayer	Yes	103022-100
XF HS Mini or XFp	V3 PS (+PDL)	8	250	2.28	200	Poly-D-Lysine on PS	2-D, suspension and monolayer	Yes	103721-100
XF HS Mini	PS	8	250	0.71	200	PS	High sensitivity* 2-D; suspension and monolayer	Yes	103725-100
XF HS Mini	PS (+PDL)	8	250	0.71	200	Poly-D-Lysine on PS	High sensitivity* 2-D; suspension and monolayer	Yes	103727-100
XF Pro	V3 PS	96	250	2.28	200	PS	2-D, monolayer	Yes	103774-100
XF96 or XF Pro	V3 PS (M Type plate)	96	250	2.28	200	PS	2-D, monolayer	Yes	103794-100
XF96 or XF Pro	V3 PS (+PDL)	96	250	2.28	200	Poly-D-Lysine on PS	2-D, suspension and monolayer	Yes	103799-100


PS = Polystyrene

* = XF HS Miniplate enables fewer cells per well.


Specifications for Cell Culture Microplates

XFe96 Cell Culture Microplate


Dimensions are [mm] in.

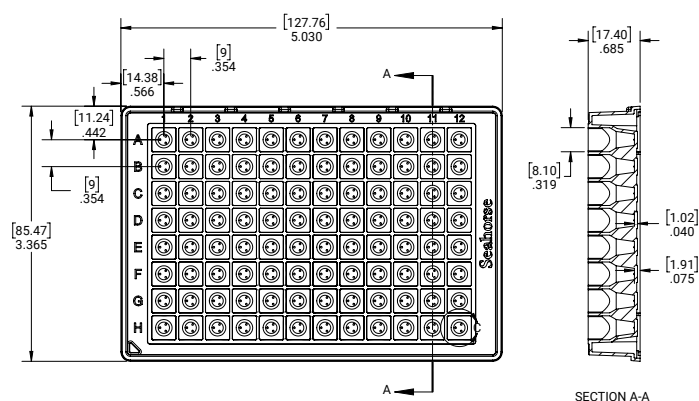


XFp Cell Culture Miniplate

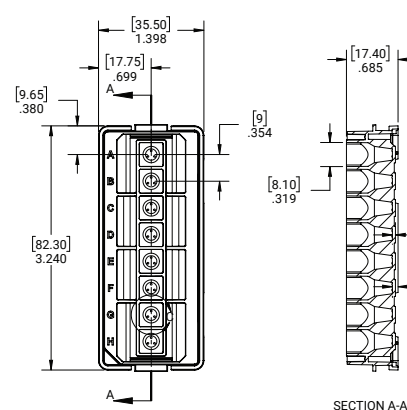

XFe24 Cell Culture Microplate

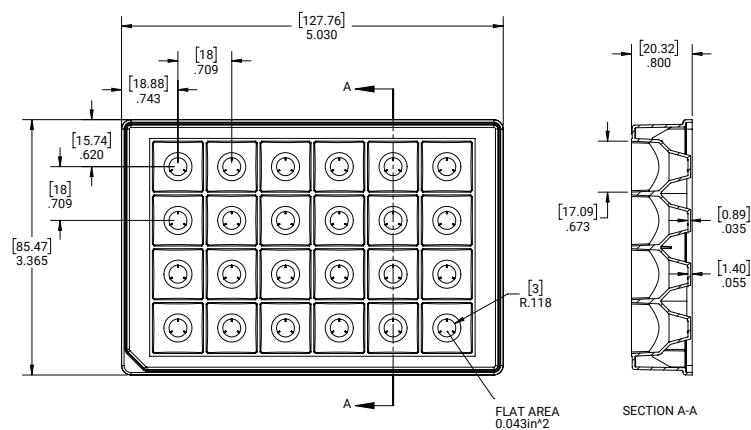

XF HS Miniplate

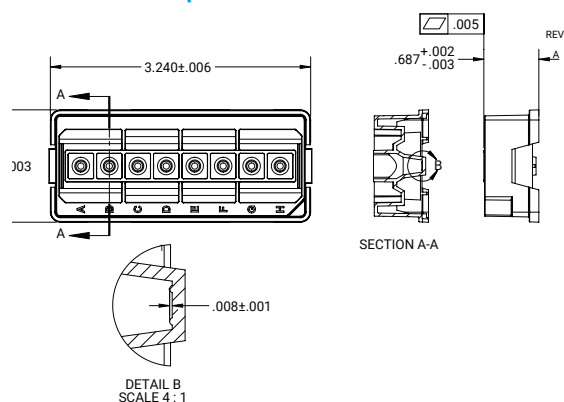

All plate dimensions can be found online
www.agilent.com/chem/seahorse-xf-plate-dimensions

Cell culture microplate dimensions

Plate Type	Max Volume (μL)	Well Depth (mm)	Well Diameter Top/Bottom (mm)	Plate Length (mm)	Plate Width (mm)	Plate Height (mm)	A1 Row Offset (mm)	A1 Column Offset (mm)	Well-to-Well Row Center Spacing (mm)	Well-to-Well Column Center Spacing (mm)	Well Bottom Elevation (mm)	Well Bottom Thickness (mm)	Well Bottom Area (mm²)
XFe96	250	15.49	8.1/3.81	127.76	85.47	17.4	11.24	14.38	9	9	0.89	1.02	11.40
XFe24	1000	18.92	17.09/6	127.76	85.47	20.32	15.74	18.88	18	18	0.51	0.89	28.26
XFp	250	15.49	8.1/3.81	35.5	82.3	17.4	9.65	17.75	9	N/A	0.89	1.02	11.40
XFe24 Islet	1000	19.18	17.09/3.15	127.76	85.47	20.32	15.74	18.88	18	18	0.5	0.65	7.79
XFe96 Spheroid	250	15.75	8.1/3.18	127.76	85.47	17.4	11.24	14.38	9	9	0.89	0.76	7.94
XF Pro M	250	15.49	8.1/3.81	127.76	85.47	17.4	11.24	14.38	9	9	0.89	1.02	11.40
XFe96/XF Pro	250	15.49	8.1/3.81	127.76	85.47	17.4	11.24	14.38	9	9	0.89	1.02	11.40

Ordering information for Agilent Seahorse XF plastic ware

Product Description	Product Content	Part Number
XF Pro Analyzer		
Seahorse XF Pro M FluxPak	18 96-well sensor cartridges, 18 XF Pro M microplates	103775-100
Seahorse XF Pro M FluxPak Mini	6 96-well sensor cartridges, 6 XF Pro M microplates	103777-100
Seahorse XF Pro M Microplates	6 XF Pro M microplates	103774-100
Seahorse XFe96/XF Pro FluxPak	18 96-well sensor cartridges, 18 XFe96/XF Pro microplates	103792-100
Seahorse XFe96/XF Pro FluxPak Mini	6 96-well sensor cartridges, 6 XFe96/XF Pro microplates	103793-100
Seahorse XFe96/XF Pro Microplates	6 XFe96/XF Pro microplates	103794-100
Seahorse XFe96/XF Pro PDL FluxPak Mini	6 96-well sensor cartridges, 6 XFe96/XF Pro PDL microplates	103798-100
Seahorse XFe96/XF Pro PDL Microplates	6 XFe96/XF Pro PDL microplates	103799-100
Seahorse XFe96 Spheroid FluxPak	6 XFe96 sensor cartridges, 6 XFe96 spheroid culture micoplates	102905-100
Seahorse XFe96 Spheroid Plate	1 polystyrene plate, 1 mm (divot diameter)	102959-100
Seahorse XFe96 Spheroid Microplates	6 polystyrene plates, 1 mm (divot diameter)	102978-100
XFe96 Analyzers		
Seahorse XFe96/XF Pro FluxPak	18 96-well sensor cartridges, 18 XFe96/XF Pro microplates	103792-100
Seahorse XFe96/XF Pro FluxPak Mini	6 96-well sensor cartridges, 6 XFe96/XF Pro microplates	103793-100
Seahorse XFe96/XF Pro Microplates	6 XFe96/XF Pro microplates	103794-100
Seahorse XFe96/XF Pro PDL FluxPak Mini	6 96-well sensor cartridges, 6 XFe96/XF Pro PDL microplates	103798-100
Seahorse XFe96/XF Pro PDL Microplates	6 XFe96/XF Pro PDL microplates	103799-100
Seahorse XFe96 FluxPak	18 XFe96 sensor cartridges, 20 XF96 V3 PS tissue culture microplates	102416-100
Seahorse XFe96 FluxPak Mini	6 XFe96 sensor cartridges, 10 XF96 V3 PS tissue culture microplates	102601-100
Seahorse XFe96 Spheroid FluxPak	6 XFe96 sensor cartridges, 6 XFe96 spheroid culture micoplates	102905-100
Seahorse XF96 V3 PS Culture Microplates	10 polystyrene plates	101085-004
Seahorse XFe96 Spheroid Plate	1 polystyrene plate, 1 mm (divot diameter)	102959-100
Seahorse XFe96 Spheroid Microplates	6 polystyrene plates, 1 mm (divot diameter)	102978-100
Seahorse XFe96 FluxPak Mini (PDL Microplates)	6 XFe96 sensor cartridges, 6 XFe96 PDL microplates	103729-100
Seahorse XFe96 PDL Microplates	6 XFe96 PDL microplates	103730-100
XFe24 Analyzers		
Seahorse XFe24 FluxPak	18 XFe24 sensor cartridges, 20 XF24 V7 PS tissue culture microplatess	102340-100
Seahorse XFe24 FluxPak Mini	6 XFe24 sensor cartridgess, 10 XF24 V7 PS tissue culture microplates	102342-100
Seahorse XFe24 Islet Capture FluxPak	6 XFe24 sensor cartridges, 6 XF24 islet capture culture microplates	103518-100
Seahorse XF24 V7 PS Culture Microplates	10 polystyrene plates	100777-004
Seahorse XF24 V28 PS Culture Microplates	10 polystyrene plates	100882-004
Seahorse XF24 Islet Capture Microplates	6 polystyrene plates	101122-100
XFp Analyzers		
Seahorse XFp FluxPak	12 XFp sensor cartridges, 12 XFp PS tissue culture microplates	103022-100
Seahorse XFp Cell Culture Miniplates	12 polystyrene plates	103025-100
Seahorse XFp FluxPaks (PDL Miniplates)	12 8-well sensor cartridges, 12 XFp miniplates	103721-100
Seahorse XFp PDL Miniplates	12 XFp PDL miniplates	103722-100
XF HS Mini Analyzers		
Seahorse XFp FluxPak	12 XFp PS tissue culture microplates	103022-100
Seahorse XFp Cell Culture Miniplates	12 polystyrene plates	103025-100
Seahorse XFp FluxPak (PDL Miniplates)	12 8-well sensor cartridges, 12 XFp PDL miniplates	103721-100
Seahorse XFp PDL Miniplates	12 XFp PDL miniplates	103722-100
Seahorse XF HS Mini FluxPak	12 8-well sensor cartridges, 12 XF HS miniplates	103723-100
Seahorse XF HS Mini FluxPak (PDL Miniplates)	12 8-well sensor cartridges, 12 XF HS PDL miniplates	103724-100
Seahorse XF HS Miniplates	12 XF HS miniplates	103725-100
Seahorse XF HS PDL Miniplates	12 8-well sensor cartridges, 12 XF HS PDL miniplates	103727-100

PS = Polystyrene

PDL = Poly-D-Lysine

Learn more about Agilent Seahorse XF technology:

www.agilent.com/chem/discoverXF

Find a local Agilent customer center:

www.agilent.com/chem/contactus

U.S. and Canada

1-800-227-9770; select option 3, then 8

cellanalysis.support@agilent.com

Europe

UK: **0800 096 7632**

Germany: **0800 180 66 78**

Netherlands: **0800 022 7243**

Other EU: **+45 3136 9878**

cellanalysis.support@agilent.com

Asia Pacific

cellanalysis.support@agilent.com

For Research Use Only. Not for use in diagnostic procedures.

RA44491.5509027778

This information is subject to change without notice.

© Agilent Technologies, Inc. 2021, 2022
Published in the USA, January 26, 2022
5991-8657EN

