

Authors

Jean Daniel Berset Water and Soil Protection Laboratory (WSPL), Bern, Switzerland

Emma Rennie Agilent Technologies, Inc., Santa Clara, CA, USA

Thomas Glauner Agilent Technologies GmbH, Waldbronn, Germany

Screening and Identification of Emerging Contaminants in Wastewater Treatment Plant Effluents Using UHPLC/Q-TOF MS and an Accurate Mass Database and Library

Application Note

Environmental, Emerging contaminants, Water, Accurate mass screening, pesticides, PPCPs

Abstract

This application note describes the creation of an accurate mass library for relevant environmental contaminants and its application for the screening of pesticides, pharmaceuticals and personal care products, their metabolites, and transformation products in effluents of municipal waste water treatment plants (WWTPs). An Agilent 1290 Infinity LC System coupled to an Agilent 6550 iFunnel Q-TOF LC/MS System was operated in positive and negative electrospray mode using Dual Spray Agilent Jet Stream Technology. Accurate mass spectra were acquired for a large collection of potential environmental contaminants in either one or both ionization modes, and for all relevant ion species.

Four WWTPs in central Europe were sampled over 3.5 months, and samples were analyzed using a Target Screening approach including reference standards for 390 contaminants. In addition, an extended suspect screening was done using all compounds included in the Agilent MassHunter Water Screening Personal Compound Database and Library (PCDL) as the suspect list. Samples were measured using All lons MS/MS as well as auto MS/MS acquisition with an inclusion list. The results clearly show the value of both acquisition modes combined with an efficient data analysis workflow, as well as the inherent value of the Water Screening PCDL for the surveillance screening of potential contaminants in complex environmental samples.



Introduction

Environmental regulations throughout the world currently focus on monitoring a limited number of well-known compounds that are assumed to be responsible for significant ecological and human health related risks [1]. As a consequence, there is a tendency to further reduce the maximum allowable concentrations of the environmental quality standards (EQS) for those priority pollutants. In Europe, for example, EU directive 2013/39/EU specifies EQS for 41 priority substances or substance groups amongst which are pesticides, perfluorocarbons (PFCs), nonyl- and octylphenol, polybrominated flame retardants, dioxins and PCBs, polyaromatic hydrocarbons (PAHs), and heavy metals [2]. However, these priority pollutants represent only a small fraction of the anthropogenic chemicals that are used, and consequently found in the environment. Apart from the legal requirement for surveillance screening for potential contaminants in surface waters in the EU, there is a growing interest to collect occurrence data for contaminants of emerging concern. This is reflected in a number of country-specific government regulations, such as the EPA method 1694, on the analysis of pharmaceuticals and personal care products in water, soil, and sediments [3], and in the US EPA Contaminant Candidate List (CCL) fourth edition from 2015 [4].

Most anthropogenic contaminants enter the aquatic environment as a result of incomplete removal in wastewater treatment plants (WWTPs). At the same time, degradation and transformation products are formed during biological wastewater treatment. These products are typically not monitored, and often are not even known [5,6]. Surface waters downstream of WWTPs contain innumerable anthropogenic contaminants and their transformation products at trace levels. When river water is subsequently used as the source for drinking water production, referred to as unplanned indirect potable reuse, comprehensive data on the chemical quality of the water are essential. This becomes even more important as planned indirect and direct potable reuse becomes a common practice [7] due to climate change, population growth, and water scarcity.

Targeted analytical methods are increasingly complemented by untargeted acquisition methods using high resolution accurate mass Q-TOF LC/MS due to comprehensive screening requirements in current environmental regulations as well as increasing interest in the occurrence of contaminants of emerging concern. Agilent Q-TOF LC/MS instruments allow for full spectrum acquisition with high sensitivity and at a very high data rate. This is essential to obtain information on molecular ions, isotope patterns, and fragments with a single large volume direct injection of water samples. For most of the LC/MS amenable compounds specified in the EU Water Framework Directive, the EU Drinking Water Directive, or the EPA Clean Water Act, method detection limits in the low ng/L range can be achieved with this type of Q-TOF LC/MS method [8].

Three different approaches towards substance identification are used in the environmental context in combination with accurate mass LC/MS: Target, Suspect, and Non-target Screening.

In Target screening, a reference standard is measured with the same analytical method and within the same worklist as the unknown sample, so that retention time, accurate mass, and fragment information can be directly compared, and (semi-)quantitative results are obtained.

Suspect screening describes a workflow where compound database and MS/MS library information of expected contaminants are used to screen for the suspected substances. While no reference standard is available, identification of compounds with increasing confidence is obtained by comparing accurate mass and isotope patterns, retention times, and coeluting fragments, or by a spectral MS/MS library comparison.

Non-target screening aims at the identification of all remaining components detected in a sample where no structural information is available. Often, non-target screening includes the statistical comparison of two or more samples or sample groups for data reduction. A full identification considers the accurate mass, isotope pattern, and MS/MS spectrum information, and includes spectral library comparison for known compounds, and spectral similarities and in-silico fragmentation tools such as Agilent MassHunter Molecular Structure Correlator (MSC) for unknowns.

Compound databases containing accurate mass MS/MS spectral data for all expected contaminants, as well as those theoretically predicted, and their transformation products have been proven to assist in the identification of potentially relevant compounds and support efficient data analysis in all three screening workflows. The Agilent MassHunter Water Screening Personal Compound Database and Library (PCDL) contains a relevant list of more than 1,400 environmental contaminants including pesticides, pharmaceuticals, personal care products, large volume industrial chemicals, and their transformation products. This PCDL includes all compounds currently regulated in the US, EU, and China, and further compounds that have either been previously detected in the environment or are likely to be detected due to their production amount or widespread use. Searchable fields containing compound information including compound class and regulation tags make it easy to create subsets of the Water Screening PCDL for target and suspect screening. The Water Screening PCDL contains more than 1,000 compounds with accurate mass MS/MS spectra, which directly supports the Agilent All lons MS/MS workflow for Target and Suspect screening, and greatly increases confidence in the identification of potential contaminants [9].

This application note describes the screening and (semi-)quantification of contaminants in WWTP effluents using an Agilent 1290 Infinity UHPLC coupled to the highly sensitive Agilent 6550 iFunnel Q-TOF LC/MS system. Four WWTPs located in agricultural and rural areas in Switzerland were sampled over 3.5 months, covering the main pesticide application period from March to late June. Comprehensive screening of potential water contaminants revealed the chemical characteristics of the different treatment and different catchment areas. A subset of the Water Screening PCDL, containing roughly 390 compounds, was used for Target screening and (semi-)guantification. For Suspect screening, the entire comprehensive Water Screening PCDL was used to find and identify compounds with high confidence. Examples of the two different screening strategies used are shown. First, the Q-TOF was operated in the All lons MS/MS mode with three collision energies. The All lons technique features an easy setup of the acquisition method with verification of the potential contaminants using the MS/MS spectral library to produce chromatographic coelution of the precursor and product ions. An additional approach was used for compounds for which either no library spectrum was available, or which were not verified due to a low coelution score. Their precursor masses were added to an inclusion list for an Auto MS/MS method using a second injection. The obtained spectra were compared to the MS/MS library for compound identification. This workflow can also be combined with a discovery run operated in TOF mode to obtain the target list of suspects.

Experimental

Reagents and standards

All reagents and solvents were HPLC or LC/MS grade. Acetonitrile, methanol, and acetic acid were purchased from Fluka (Sigma-Aldrich, Buchs, Switzerland). Ammonium acetate was purchased from VWR International (Darmstadt, Germany). Ultrapure water was produced using a Milli-Q Integral system equipped with a 0.22-µm point-of-use membrane filter cartridge (EMD Millipore, Billerica, MA, USA). For acquisition of accurate mass MS/MS spectra, most individual reference compounds were purchased from Sigma-Aldrich (Buchs, Switzerland) or from VWR International (Darmstadt, Germany). Reference standards for transformation products and residues were purchased from Ehrenstorfer (LGC Standards, Wesel, Germany). Estimation of retention times was performed using mixed standard solutions of pesticides, pharmaceuticals, and drugs of abuse, which were part of the LC/MS Pesticide Comprehensive Test Mix (p/n 5190-0551), the LC/MS Forensic Toxicology Comprehensive Test Mix (p/n 5190-0555), or were provided by research groups.

Stock standard solutions were prepared by dissolving the reference compounds in isopropanol, acetonitrile, methanol, water, or mixtures thereof, depending on the physicochemical properties of the substance. For calibration and spiking experiments, the comprehensive test mixtures for pesticides, pharmaceuticals, and drugs of abuse were combined to a multi-analyte working solution. Stock standard solutions as well as the multi-analyte working solution were stored until use at -20 °C. Calibration samples were prepared by dilution of the working solution with tap water.

Sample preparation

Effluents of four different wastewater treatment plants in central Europe were collected as 14-day composite samples over 3.5 months (March to June). The catchment areas of three of the WWTPs are agriculturally dominated (AG, AI, AL), whereas the other one is located in an urban area (AZ). One of the WWTPs also receives some industrial wastewater (AL). Efficient nitrification-denitrification is observed in three of the WWTPs (AG, AL, AZ) and, thus, a better elimination of trace contaminants can be expected.

Samples were filtered using glass fiber filters, and were stored at -20 °C. Immediately before measurement, samples were thawed, and an aliquot was transferred to an HPLC vial.

LC/Q-TOF MS analyses

Separation was carried out using an Agilent 1290 Infinity UHPLC system consisting of:

- Agilent 1290 Infinity Binary Pump (G4220A)
- Agilent 1290 Infinity High Performance Autosampler (G4226A) equipped with a large volume injection kit (G4216-68711)
- Agilent 1290 Infinity Thermostatted Column compartment (G1316C)

The UHPLC system was coupled to an Agilent G6550A iFunnel Quadrupole Time-of-Flight LC/MS System equipped with a Dual Spray Agilent Jet Stream electrospray ionization source. The Q-TOF LC/MS instrument was operated with:

- Agilent MassHunter Acquisition rev. B.06.01 in 2 GHz extended dynamic range mode with positive or negative electrospray ionization (ESI) with two different methods:
 - All lons MS/MS acquisition with 3 scans/sec with two discrete collision energies
 - Auto MS/MS acquisition with a data rate of 5 scans/sec in MS and 5 scans/sec in MS/MS

The use of two collision energies in the All lons acquisition resulted in alternating spectra with a low energy channel containing the precursor ion, and two high energy channels containing the precursor and product ions.

Reference mass ions were delivered using an Agilent Infinity 1260 Isocratic pump (G1310B) operated with a flow rate of 1.0 mL/min, and using a 1-in-100 flow splitter (G1607-60000). The final flow rate going to the reference sprayer was 10 μ L/min.

Table 1 shows the chromatographic conditions, and Table 2 shows the major MS conditions.

Data were evaluated using MassHunter Qualitative Analysis software B.07.00. Positive identifications of water contaminants were reported if the compound was detected in the accurate mass MS data by the Find-by-Formula data mining algorithm with a mass error below 5 ppm, and with a sufficient score (including isotope abundance and isotope spacing). A retention time window of ±1 minute was specified for peak detection to compensate for retention time shifts due to matrix variability.

Table 1. Chromato	graphic Conditions					
Parameter	Value					
UHPLC column	Agilent ZORBAX RRHD SB-Aq, 2.1 × 150 mm, 1.8 μm (p/n 859700-914)					
Column temperature	40 °C	40 °C				
Mobile phase	A) 1 mM NH ₄ acetate + 0.1% acetic acid B) 0.1% acetic acid in acetonitrile					
Gradient program	Time (min) 0.0 2.0 14.0 16.0 19.0 19.5	% B 0 0 98 98 0 0				
Stop time	20.0 minutes					
Post time	3.0 minutes					
Flow rate	0.40 mL/min					
Injection volume	100 µL					

Table 2. Major MS Conditions

Parameter	Value				
Gas temperature	160 °C				
Gas flow	16 L/min				
Nebulizer	30 psig				
Sheath gas temperature	350 °C				
Sheath gas flow	12 L/min				
	Positive	Negative			
Capillary voltage	4,500 V	3,500 V			
Nozzle voltage	500 V	1,000 V			
Reference mass correction	121.05087 922.00980	112.98559 1,033.98811			
All ions MS/MS					
Mass range	50 to 1,200 amu				
Scan rate	3 spectra/s				
Collision energies	0-20-40 V				
Auto MS/MS					
MS Mass range	100 to 1,200 amu				
MS/MS Mass range	50 to 1,200 amu				
Scan rate	5 spectra/s 5 spectra/s	(MS) (MS/MS)			
Collision energy	20 V				

Creation of the Agilent MassHunter Water Screening PCDL

Accurate mass spectra of single-analyte solutions were acquired with flow injection or using a short column in targeted MS/MS mode with collision energies of 10, 20, and 40 V. If precursor ion stability required higher collision energies, additional spectra were acquired in a second run. Typically, MS/MS spectra were acquired for the [M+H]⁺ and [M-H]- ion species for each analyte. If highly abundant additional adduct ion species were observed, accurate mass MS/MS spectra were also acquired for the [M+NH₄]⁺, [M+Na]⁺, or [M+Cl]⁻ species. In either positive or negative ionization mode, meaningful MS/MS spectra were acquired for more than 1,000 relevant water contaminants. For many compounds, MS/MS library spectra were captured in both ionization modes, and for more than one precursor ion species. To eliminate mass assignment errors, fragment masses in the acquired spectra were compared to the theoretical fragment formulas, and all ion peaks were

corrected to their theoretical masses. All MS/MS spectra were curated for spectral noise, and a minimum base peak threshold was applied to ensure good ion statistics for all fragment ions. The corrected spectra were included in the Agilent Water-Screening Personal Compound Database and Library (G6882CA).

The Water Screening PCDL was then used for the screening and identification of environmental contaminants in the effluents of four WWTPs. Moreover, by analyzing mixed standard solutions of pesticides, pharmaceuticals, and drugs of abuse with the given UHPLC method, retention time information was added to 390 compounds, thus adding in retention time to the identification score and increasing identification confidence.

Figure 1 shows a screen capture of the MassHunter PCDL Manager software along with the accurate mass MS/MS spectrum of the antiviral drug aciclovir acquired in negative ionization mode with a collision energy of 20 V.

	🔛 MassHunter PCDL Manager for Forensics and Toxicology - C:\MassHunter\PCDL\WaterScreening_AM_PCDL.cdb												
÷F	ile Edit View PCDL Links Help												
	Find Spectra 🎒 📃 🗋 📴 😹 🛗 🥹												
	Single Sarah Patah Sarah	Patch Summan	Edit Como	unda	Con	ustral Casurals	Browse	Spectra	Edit Spacetra				
	Single Search	bateri Summary	Luit Comp	Julius	Spe	Scial Search			Graphin Mass List				
ſ	Mass	lon polarity:	(Any)						Chapme Mass List				
	Precursor ion:		(***)						□ Library spectrum □ 110 1 150 04213				
	Tolerance: 200 💿 ppm 💿 mDa	lonization mode:	(Any)		•				5 100 - 100.000 - 224 0789				
1	Collision energy												
									00				
	Tolerance: 2.0 eV								20				
									70-				
S	pectra for compound: Aciclovir (Acyclovir)								60-				
	Compound Name	Ion Species	Precursor lo	CE (V)	Pol	arity Ioniz	ation Instrum	nent	50-92.02542				
	Aciclovir (Acyclovir)	(M-H)-	224.078	1	10 Neg	ative ESI	QTOF		40-				
Þ	Aciclovir (Acyclovir)	(M-H)-	224.078	1	20 Neg	ative ESI	QTOF		30-				
	Aciclovir (Acyclovir)	(M-H)-	224.078	1	40 Neg	ative ESI	QTOF		20- 66.00977 19.58				
	Aciclovir (Acyclovir)	(M+H)+	226.093	7	10 Posi	tive ESI	QTOF		10-				
	Aciclovir (Acyclovir)	(M+H)+	226.093	7	20 Posi	tive ESI	QTOF						
	Aciclovir (Acyclovir)	(M+H)+	226.093	7	40 Posi	tive ESI	QTOF		20 40 60 80 100 120 140 160 180 200 220 2	240			
									m/z				
	Si	ngle Search Res	ults: 1193	hits									
	Compound Name	Fomula	Mass	Anion	Cation	RT (min)	CAS	ChemSpider	IUPAC Name Spectra	^			
	Paraxanthine	C7H8N4O2	180.06473				611-59-6	4525	1,7-Dimethyl-3,7-dihydro-1H-purine-2,6-dione 6				
	Saccharin	C7H5NO3S	182.99901			6.755	81-07-2	4959	1.2-Benzothiazol-3(2H)-one 1.1-dioxide 6				
	Acephate	C4H10NO3PS	183.01190				30560-19-1	1905	O,S-Dimethyl acetylphosphoramidothioate 6				
	Sulfacetamide	C8H10N2O3S	214.04121				144-80-9	<u>5129</u>	N-[(4-Aminophenyl]sulfonyl]acetamide 6				
	Primidone	C12H14N2O2	218.10553			7.407	125-33-7	4740	5-Ethyl-5-phenyldihydro-4,6(1H,5H)-pyrimidinedione 6				
	Metaxalone	C12H15NO3	221.10519				1665-48-1	14709	5-[(3,5-Dimethylphenoxy)methyl]-1,3-oxazolidin-2 6				
	Tapentadol	C14H23NO	221.17796				175591-23-8	8013742	3-[(1R,2R)-3-(dimethylamino)-1-ethyl-2-methylprop 6				
	Acetazolamide	C4H6N4O3S2	221.98813				<u>59-66-5</u>	<u>1909</u>	N-(5-Sulfamoyl-1,3,4-thiadiazol-2-yl)acetamide 6				
►	Aciclovir (Acyclovir)	C8H11N5O3	225.08619				<u>59277-89-3</u>	<u>1945</u>	2-Amino-9-[(2-hydroxyethoxy)methyl]-1,9-dihydro-6 6				
	Naproxen	C14H14O3	230.09429			11.237	22204-53-1	<u>137720</u>	(2S)-2-(6-Methoxy-2-naphthyl)propanoic acid 6				
	Diuron	C9H10Cl2N2O	232.01702			10.810	<u>330-54-1</u>	3008	3-(3,4-Dichlorophenyl)-1,1-dimethylurea 6				
	Nalidixic acid	C12H12N2O3	232.08479				<u>389-08-2</u>	<u>4268</u>	1-Ethyl-7-methyl-4-oxo-1,4-dihydro-1,8-naphthyridi 6				
	Salbutamol (Albuterol)	C13H21NO3	239.15214				18559-94-9	<u>1999</u>	2-(Hydroxymethyl)-4-{1-hydroxy-2-{(2-methyl-2-pro 6				
	Mefenamic acid	C15H15NO2	241.11028			12.855	<u>61-68-7</u>	<u>3904</u>	2-[(2,3-Dimethylphenyl)amino]benzoic acid 6				
	Cimetidine	C10H16N6S	252.11572				<u>51481-61-9</u>	2654	1-Cyano-2-methyl-3-(2-{[(4-methyl-1H-imidazol-5-yl) 6				
	Fenbufen	C16H14O3	254.09429	(T)			36330-85-5	3218	4-(4-BiohenvlvI)-4-oxobutanoic acid 6	-			

Figure 1. Agilent MassHunter PCDL Manager software showing the Agilent MassHunter Water Screening PCDL and the accurate mass MS/MS spectrum of aciclovir acquired in negative ionization mode with a collision energy of 20 V.

Results and Discussion

Target screening for pesticides, pharmaceuticals, and personal care products

For the target screening and (semi-)quantification of pesticides, pharmaceuticals, and personal care products in the WWTP effluents, samples were measured in positive and negative All lons MS/MS mode with collision energies of 0, 20, and 40 V. A subset PCDL with 390 entries was created from the Water Screening PCDL including all compounds for which reference standards were available. Data for the calibration samples were initially evaluated in the MassHunter Qualitative Analysis Software (B.07.00) using the Find-by-Formula data mining algorithm with a mass error of ± 5 ppm and a retention time window of ± 0.5 minutes compared to the expected retention time. In positive mode, [M+H]⁺ and [M+NH,]⁺ species, and in negative mode, [M-H]⁻ species have been considered as charge carriers since they are represented by the majority of accurate mass MS/MS spectra included in the PCDL. Identification was done by Fragment Confirmation using the Water Screening PCDL as the fragment ion source, and evaluating the five most specific ions from the MS/MS spectral library. In the All lons

MS/MS software workflow, fragment ions that show perfect coelution with the molecular ions were identified and the ideal collision energies and relative ratios were detected. This information was passed to the MassHunter Quantitative Analysis software for (semi-)quantification and batch review using accurate mass for molecular ions and fragments as well as isotope pattern matching for compound identification with high confidence.

From the 390 targeted compounds, 315 were detected in positive mode, and 75 were detected in negative mode. With a direct injection of 100 μ L of water into the UHPLC Q-TOF MS system, more than 60% of the compounds could be quantified at or below 10 ng/L in the spiked tap water samples. Another 35% of the compounds were detected between 10 and 100 ng/L, and just 5% of the compounds were only detected at concentrations of 200 ng/L or higher. For most targeted compounds, one or more specific fragment ions could be used as qualifier ions, and generally mass accuracy of the molecular ions and fragments was better than 5 ppm. Figure 2 shows the extracted ion chromatograms, MS peak spectra and calibration curves for some examples from the lists of priority pollutants (EU) and US EPA method 1694.



Figure 2. EIC chromatograms of molecular ion and fragments, MS peak spectra and calibration curves for erythromycin (positive mode), ibuprofen (negative mode), diclofenac (negative mode), and quinoxyfen (positive mode).

Applying this target screening method to the samples from the WWTP effluents allowed the (semi-)quantification of a large variety of environmental contaminants in a concentration range of a few ng/L up to several μ g/L. Pharmaceutical residues were found to dominate the effluents of the larger WWTPs (AG, AL, AZ), which all receive wastewater from about 50,000 population equivalents, while Al receives wastewater from 7,100 population equivalents. In all wastewater treatment plant effluents, X-ray contrast media were identified with estimated concentrations of up to 7 μ g/L for iomeprol, and up to 2 μ g/L for iopromide. Other common pharmaceuticals were amisulprid (up to 500 ng/L), atenolol (up to 1.7 μ g/L), metoprolol (up to 470 ng/L), and tramadol (up to $2 \mu g/L$) as well as carbamazepine, diclofenac, ibuprofen, naproxen, and sulfamethoxazole. Metabolites were identified for carbamazepine, sulfamethoxazole, and metamizole. In total, 33 pharmaceuticals and metabolites of pharmaceuticals were identified in the WWTP effluents. As expected, the largest number of these compounds were found in the effluent of AZ, which has an urban catchment area. However, concentrations were highest in effluents of Al, probably due to the limited elimination efficiency of the simple treatment.

In the effluents of all WWTPs, a total of 46 pesticides were detected. The insecticide diethyltoluamide (DEET) was found in all WWTPs with a concentration of 14 to 770 ng/L. Also present in samples of all WWTPs were the herbicides metolachlor (up to $1.1 \ \mu g/L$) and isoproturon (up to $450 \ ng/L$). The largest number of pesticides were found in the effluent of Al, which has an agriculturally-dominated catchment area. Major crops grown there include cereals, vegetables, corn, beetroot, and potatoes. The pesticides found most often, and with the highest concentrations in the effluent of this WWTP were azoxystrobin, flufenacet, linuron, metamitron, methomyl, metribuzin, propamocarb, spiroxamin, and terbuthylazine. All of these pesticides are mainly used for the above listed crops.

Figure 3 shows the normalized chromatograms for the antiepileptic drug carbamazepine and the pesticide azoxystrobin in the effluent samples from WWTP AI, spanning a time period from March to end of June. While the concentration of the pharmaceutical is more or less constant over the whole sampling period, the concentration of the pesticide increased over the growing season. This example demonstrates the different contaminant profiles for compounds that are continuously introduced, compared to those that enter the water cycle within a specific period of time.



Figure 3. Normalized EIC chromatograms of molecular ions and fragments for the antiepileptic drug carbamazepine and the pesticide azoxystrobin over the course of the sampling period (March to end of June).

Extended Suspect screening using the comprehensive Agilent MassHunter Water Screening PCDL

Based on the results from the target screening, the samples from the WWTPs AI and AZ were selected for an extended Suspect screening looking for all remaining compounds included in the Water Screening PCDL. The availability of accurate mass MS/MS information is key for the identification of potential candidates, and is either used in the All lons MS/MS workflow for the extraction and alignment of EICs of the molecular ion and characteristic fragments, or for the library matching of an acquired accurate mass MS/MS spectrum against the reference spectra in the PCDL. Both workflows are shown in Figure 4 for some of the examples identified in the WWTP effluents.

All lons MS/MS screening workflow

In the All lons MS/MS workflow, accurate mass data is collected without fragmentation in a low energy channel.

Virtually at the same time, using two high energy channels, compounds are fragmented with two different collision energies without precursor selection in very fast sequential steps, and accurate mass fragment data are recorded for both channels. When the data is analyzed using the Find-by-Formula algorithm, the Water Screening PCDL provides the precursor formula information, and compound chromatograms are extracted for all specified ion species. For putative identifications, chromatograms are automatically extracted from the high energy channels for a specified number of the most abundant fragments from the MS/MS spectra stored in the PCDL. As an example, Figure 4A shows the accurate mass library spectrum of valsartan from the Water Screening PCDL, compared to the cleaned high energy spectrum from an effluent sample from the WWTP AI (Figure 4B). The red triangles indicate the fragment ions that have been selected from the library spectrum for evaluation.



Figure 4. Accurate mass library spectrum for valsartan at a collision energy of 20 V (A) in comparison to the acquired high energy spectrum (B) from an effluent sample of the wastewater treatment plant Al (cleaned spectrum). The red triangles in the library spectrum (A) indicate automatically selected ions for the All lons MS/MS evaluation.

While the library spectrum is based on a collision energy of 20 V, the cleaned high energy spectrum combines information from both high energy channels acquired with 20 and 40 V.

By overlaying chromatograms for both precursor and fragment ions, and the calculation of a coelution score, the identity of the angiotensin receptor blocker (ARB) valsartan was confirmed. The coelution score takes into account factors such as abundance, peak shape (symmetry), peak width, and retention time. The normalized intensity ratios are plotted and made available to the user for inspection in a coelution plot. Figure 5A shows the overlay of the molecular ion chromatogram with the fragment chromatograms from the high energy channels. All five fragment chromatograms showed coelution with the precursor ion. This is also demonstrated in the coelution plot in Figure 5B. The detailed compound identification results, including the coelution scores, are shown in the compound table in Figure 5C.

Identification with high confidence is achieved when the EICs of the molecular ion and at least one or two fragments show perfect coelution, which is expressed by a coelution score of > 90 (out of 100), and the mass accuracy for the peak spectrum for both molecular ions and fragments is better than 5 ppm. Based on this rule, additional pharmaceuticals were identified in the effluent of WWTP AI, namely the ARBs candesartan, irbesartan, and losartan.



Figure 5. Overlay of precursor and fragment ion traces for valsartan in an effluent sample of the wastewater treatment plant AI (A), coelution plot (B) and compound identification results including the coelution score (C).

Other pharmaceuticals and personal care products that were not targeted in the previous workflow (target screening using a reference standard), including the compounds clarithromycin, fexofenadine, sitagliptin, celiprolol, and phenylbenzimidazole sulfonic acid were also identified. Moreover, nine further pesticides (napropamid, pyrimethanil, fenamidone, lenacil, dimethenamid, boscalid, dinoseb, fludioxonil, and penconazole) as well as perfluorooctanoic acid (PFOA) and several organophosphates (triethyl phosphate, tris(2-chloroethyl)phosphate, tributylphosphate, and triphenylphosphate) were also found. In the AZ samples, valsartan, candesartan, and irbesartan, as well as fexofenadine, clarithromycin, venlafaxine, and its metabolite desmethyl venlafaxine, citalopram, cetirizine, clopidogrel, and ritonavir were found. In addition to PFOA, perfluorononanoic acid was also found. No further pesticides were detected and confirmed in the effluent of the WWTP AZ.

Suspect screening and verification using targeted MS/MS

The verification of compounds by matching accurate mass MS/MS spectra against reference spectra from an accurate mass MS/MS library sometimes offers advantages over the All lons MS/MS workflow. This is the case if contaminants are present only at trace concentrations and, thus, signal intensities of the fragment ions are low, or if interferences occur between the low mass fragment ions and ions from the matrix. Moreover, accurate mass MS/MS library comparison after precursor isolation is considered to be the gold standard

for compound verification. In cases where there is no library spectrum available, for example, for newly identified compounds or suspected transformation products, accurate mass MS/MS spectra can be compared to theoretical fragmentation of a compound in the MassHunter Molecular Structure Correlator (MSC) software.

Typically, this workflow starts with an All Ions MS/MS acquisition discovery run, and the tentative identification of suspected compounds using the Find-by-Formula data mining algorithm with the fragment confirmation. In cases when no fragment ions can be identified or if only one unspecific fragment is observed, the compounds are selected from the MassHunter Qualitative Analysis software, and exported into a target list. In a consecutive run using the same chromatography, accurate mass MS/MS spectra for the targeted precursors are acquired. If utmost sensitivity is required, it might be beneficial to operate the Q-TOF in TOF mode for the discovery run. Using TOF mode results in a higher number of suspects and, therefore, more precursor masses for the consecutive targeted MS/MS or auto MS/MS run. However, in this instance, more of the low abundant contaminants will be detected. Data analysis for the targeted MS/MS or auto MS/MS run starts with data mining using the Find by MS/MS (target or auto) feature finding algorithm followed by library searching using the MassHunter Water Screening PCDL.

Figure 6 shows the compound chromatogram and MS1 level peak spectrum for the antidiabetic drug metformin found in the effluent of wastewater treatment plant AZ. Due to the low mass of the molecular ion and the even lower masses of the specific fragments, the compound could not be verified in the All Ions MS/MS workflow, but was confidently identified with the targeted MS/MS approach. The predominant ion species for metformin was $[M+H]^+$, and the measured m/z was in good agreement with the calculated mass (–0.7 ppm). The red boxes around the mass signals show the expected isotope ratio. The measured intensities for the monoisotopic mass signal, and the [M+1] isotope signal are in very good agreement with the theoretical pattern. However, there was an interference observed for the [M+2] signal, and therefore the target score was only 72.7 (out of 100).



Figure 6. Compound chromatogram and cleaned peak spectrum obtained by the Find-by-Formula algorithm for the antidiabetic drug metformin found in the effluent of wastewater treatment plant AZ.

The red diamond in Figure 6 indicates that MS/MS spectra have been acquired for that particular *m/z*. MS/MS spectra were extracted automatically over the peak range, and were matched against the library spectra contained in the MassHunter Water Screening PCDL. Figure 7 shows the MS/MS spectrum for metformin acquired in the wastewater treatment plant effluent (7A) in comparison to the library spectrum from the PCDL (7C). A mirror representation of the

difference spectrum is shown in panel 7B. All major fragment ions listed in the library spectrum of metformin were found in the measured spectrum within a narrow mass extraction window and in a similar ratio as in the reference spectrum for a collision energy of 20 V. Thus, the reverse search against the accurate mass library resulted in a score of 95.4 out of 100, and verified the presence of metformin in the sample. MS/MS scores were required to be above 60 for verification.



Figure 7. Comparison of the measured spectrum of metformin in the WWTP effluent sample (A) with the reference spectrum from the Agilent MassHunter Water Screening PCDL (C), and a mirror representation of the difference spectrum (B).

Figure 8 shows other compound examples that could be verified by using the targeted MS/MS workflow. Melamine was found in the effluent of AZ with a library score of 85.3. It has several industrial uses, and is formed as a metabolite of the pesticide cyromazine. Denatonium was found in the effluent of AI with a library score of 91.5. It is the most bitter compound known, and is used as bitterant in personal care products to prevent inappropriate ingestion.



Figure 8. Comparison of the measured spectrum of melamine in the WWTP effluent sample AZ (A) and of denatonium in the effluent sample AI (B) with the reference MS/MS spectra from the Agilent MassHunter Water Screening PCDL.

Conclusions

A method for the screening and (semi-)quantification of environmental contaminants in water samples was developed and applied to WWTP effluents. This method takes full advantage of the low delay volumes of the Agilent 1290 Infinity LC, and its ability to handle high backpressures in UHPLC separations to increase the chromatographic resolution. The method benefits from the sensitivity of the Agilent 6550 iFunnel Q-TOF, and from the versatile ionization capabilities of the Agilent Jet Stream ionization source.

The Agilent MassHunter Water Screening PCDL is an ideal complement for the Target and suspect screening workflows in the Agilent MassHunter Qualitative and Quantitative software. In Target Screening and (semi-)quantification, the PCDL is used to define the suite of compounds and to identify selective qualifier ions. While in suspect screening, identification with high confidence is achieved by fragment confirmation or library searching. Applying both workflows to the WWTP effluent samples revealed the characteristics of the treatment technology as well as the catchment area by the different chemical inventory of trace contaminants.

Auto MS/MS acquisition with an inclusion list and a single collision energy combined with library matching resulted in similar verification rates as the Agilent All Ions MS/MS acquisition with fragment coelution. However, it can be seen that the precursor isolation in auto MS/MS acquisition improved the identification compared to the All Ions MS/MS workflow in heavy matrix and for low mass contaminants. Conversely, All Ions MS/MS acquisition is very fast, and allows the differentiation of closely eluting isomers. Another important feature of the All Ions MS/MS workflow is that the data can be re-interrogated at a later time for compounds that were not in the scope of the analysis during measurement, allowing for retrospective data analysis for new emerging contaminants without the need to reacquire data from old samples.

References

- P. Gago-Ferrero, *et al.* Extended Suspect and Non-Target Strategies to Characterize Emerging Polar Organic Contaminants in Raw Wastewater with LC-HRMS/MS. *Environ. Sci. & Technol.* 49(20), 12333-12341 (2015).
- Directive 2013/39/EU of the European Parliament and of the Council of 12 August 2013 amending Directives 2000/60/EC and 2008/105/EC as regards priority substances in the field of water policy.
- US EPA Method 1694 (2007) Pharmaceuticals and Personal Care Products in Water, Soil, Sediment, and Biosolids by HPLC/MS/MS. http://www.epa. gov/sites/production/files/2015-10/documents/ method_1694_2007.pdf (accessed 14 December 2015).
- US EPA Contaminant Candidate List (CCL) and Regulatory Determination: Draft CCL 4 Chemical Contaminants. http://www.epa.gov/ccl/chemical-contaminants-ccl-4 (accessed 14 December 2015).
- M. J. Gómez, et al. Rapid automated screening, identification and quantification of organic micro-contaminants and their main transformation products in wastewater and river waters using liquid chromatography-quadrupole-time-of-flight mass spectrometry with an accurate-mass database. J. of Chromatog. A 1217(45), 7038-7054 (2010).
- T. Letzel, *et al.* LC–MS screening techniques for wastewater analysis and analytical data handling strategies: Sartans and their transformation products as an example. *Chemosphere* **137**, 198-206 (2015).
- C. Rodriguez, *et al.* Indirect Potable Reuse: A Sustainable Water Supply Alternative. *Int. J. Environ. Res. Public Health* 6, 1174-1209 (2009).
- D-H. D. Yang, *et al.* Sensitive Screening of Pharmaceuticals and Personal Care Products (PPCPs) in Water Using an Agilent 6545 Q-TOF LC/MS System. *Agilent Technologies Application Note*, publication number 5991-5954EN (2015).
- E. L. Schymanski, *et al.* Non-target screening with high-resolution mass spectrometry: critical review using a collaborative trial on water analysis. *Anal. Bioanal. Chem.* **407(21)**, 6237-6255 (2015).

For More Information

These data represent typical results. For more information on our products and services, visit our Web site at www.agilent.com/chem.

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2016 Printed in the USA March 29, 2016 5991-6627EN

