

PicoMaxx High Fidelity PCR Master Mix

Instruction Manual

Catalog #600650 (100 reactions) Revision C.0

For Research Use Only. Not for use in diagnostic procedures. 600650-12



LIMITED PRODUCT WARRANTY

This warranty limits our liability to replacement of this product. No other warranties of any kind, express or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Agilent. Agilent shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

ORDERING INFORMATION AND TECHNICAL SERVICES

Email

techservices@agilent.com

World Wide Web

www.genomics.agilent.com

Telephone

Location	Telephone	
United States and Canada	800 227 9770	
Austria	01 25125 6800	
Benelux	02 404 92 22	
Denmark	45 70 13 00 30	
Finland	010 802 220	
France	0810 446 446	
Germany	0800 603 1000	
Italy	800 012575	
Netherlands	020 547 2600	
Spain	901 11 68 90	
Sweden	08 506 4 8960	
Switzerland	0848 8035 60	
UK/Ireland	0845 712 5292	
All Other Countries	Please visit <u>www.agilent.com/genomics/contactus</u>	

PicoMaxx High Fidelity PCR Master Mix

MATERIALS PROVIDED

Materials provided	Quantity	
PicoMaxx 2× master mix ^a	2.5 ml (100, 50-µl reactions)	

^a The total Mg²⁺ concentration present in the final $1 \times$ dilution of the $2 \times$ PicoMaxx master mix is 2 mM. The total dNTP concentration present in the final $1 \times$ dilution is 800 μ M (200 μ M each dNTP).

Storage: The PicoMaxx $2\times$ master mix should be stored at -20° C upon receipt. After thawing, store at 4° C. Once thawed, full activity is guaranteed for 6 months.

NOTICES TO PURCHASER

Use of this product is licensed under U.S. Patent Nos. 5,587,287 and corresponding foreign patents Europe, Hong Kong, and Japan

Limited Label License: The purchase of this product conveys to the purchaser only the limited, non-transferable right to use the purchased quantity of the product for the purchaser's own research by the purchaser only under patent claims outside the U.S. corresponding to the expired U.S. Patent No. 5,436,149 (claims 6-16). No rights are granted to the purchaser to sell, modify for resale or otherwise transfer this product, either alone or as a component of another product, to any third party. Takara Bio reserves all other rights, and this product may not be used in any manner other than as provided herein. For information on obtaining a license to use this product for purposes other than research, please contact Takara Bio Inc., Seta 3-4-1, Otsu, Shiga 520-2193, Japan (Fax +81-77-543-9254).

NOTICE TO PURCHASER: LIMITED LICENSE

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 6,258,569, 6,171,785, 6,127,155, 6,030,787, 5,994,056, 5,876,930, 5,804,375, 5,789,224, 5,773,258 (claims 1 and 6 only), 5,723,591, 5,677,152 (claims 1 to 23 only), 5,618,711, 5,538,848, and claims outside the US corresponding to expired US Patent No. 5,079,352. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

INTRODUCTION

The PicoMaxx High Fidelity PCR Master mix* is a 2× formulation of the PicoMaxx high fidelity PCR enzyme blend, an optimized PCR reaction buffer, magnesium, and dNTPs. The master mix includes a blend of cloned *Taq* and *Pfu* DNA polymerases** and Agilent's ArchaeMaxx polymerase-enhancing factor. Together with the optimized buffer in the master mix, this enzyme blend provides maximum PCR sensitivity and efficiency. The PicoMaxx high fidelity PCR master mix successfully amplifies even where starting material is limited, and reliably produces high PCR product yields on a wide variety of templates up to 10 kb.

The PicoMaxx high fidelity PCR master mix is formulated with antibodies that inhibit polymerase activity until cycling begins, promoting high specificity and reducing background.

PCR PROTOCOL

1. Add the PCR reaction components in order while mixing gently. Table I provides an example reaction mixture for the amplification of a typical single-copy chromosomal target. The recipe listed is for one reaction and must be adjusted for multiple samples.

Component	Amount per reaction
Distilled water (dH_2O)	22 μl
Primer #1 (100 ng/µl)∝	1 μΙ
Primer #2 (100 ng/µl)∝	1 μl
DNA template (100 ng/µl) ^b	1 μl
PicoMaxx 2 $ imes$ master mix	25 μl
Total reaction volume	50 μl

TABLE I Reaction Mixture for a Typical PCR Amplification

^a Primer concentrations between 0.3 and 0.5 μM are recommended (this corresponds to 100–200 ng for typical 18- to 25-mer oligonucleotide primers in a 50-μl reaction volume).

^b The amount of DNA template required varies depending on the type of DNA being amplified. Generally 100–200 ng of genomic DNA template is recommended. To amplify low-complexity targets (for example, lambda DNA or cloned DNA), we recommend using 5–20 ng of template. Excess template DNA can inhibit the PCR reaction.

* U.S. Patent Nos. 6,734,293, 6,444,428, 6,183,997, and 5,948,663.

Revision C.0

^{**} U.S. Patent No. 5,948,663.

- 2. Aliquot 50 μl of the reaction mixture into the appropriate number of sterile thin-wall PCR tubes or standard 0.5-ml microcentrifuge tubes.
- 3. Perform PCR using optimized cycling conditions. Suggested cycling parameters are indicated in Table II.
- 4. Analyze the PCR amplification products on a 0.7–1.0% (w/v) agarose gel.

Segment	# of Cycles	Temperature	Duration
1	1	95°C	2 minutes
2	30–35	95°C	40 seconds
		Primer T _m – 5°C ^c	30 seconds
		72°C	1 minute per kb PCR target
3	1	72°C	10 minutes

TABLE II PCR Cycling Parameters for a Typical PCR Amplification^{*a,b*}

^a Thin-wall PCR tubes are highly recommended.

^b Optimized cycling parameters are not necessarily transferable between thermal cyclers designed by different manufacturers; therefore, each manufacturer's recommendations for optimal cycling parameters should be consulted.

^c The annealing temperature may require optimization. Typically annealing temperatures will range between 55° and 72°C.¹

TROUBLESHOOTING

Observation	Suggestion(s)
No product or low yield	Increase extension time to 90 seconds per kb of PCR target.
	Use intact and highly purified DNA templates.
	Increase the amount of full-length intact DNA template, adjust the ratio of primer versus template, and/or increase the number of cycles up to a maximum of 40 cycles to optimize yield of the desired product.
	Remove extraneous salts from the PCR primers and DNA preparations.
	For some targets, increasing the dNTP concentration by supplementing with additional dNTPs may be beneficial.
	Check the melting temperature, purity, GC content, and length of the primers.
	Use the recommended primer concentrations between 0.3 and 0.5 μ M (generally 100–250 ng for typical 20- to 30-mer oligonucleotide primers in a 50- μ l reaction volume).
	Primer pairs with matching primer melting temperatures (T _m) and complete complementarity between the primer and template are recommended.
	If using existing primer pairs, the annealing temperatures may require optimization.
	Denaturation times of 30–60 seconds at 94–95°C are usually sufficient, while longer denaturation times may damage the DNA template. Use the shortest denaturation time compatible with successful PCR on the thermal cycler.
	Use thin-wall PCR tubes. These PCR tubes are optimized to ensure ideal contact with the block to permit more efficient heat transfer and to maximize thermal-cycling performance.
Artifactual smears or multiple	Increase the annealing temperature in 4°C increments.
bands	Optimize the cycling parameters specifically for the primer-template set and the thermal cycler used.

REFERENCES

1. Innis, M. A., Gelfand, D. H., Sninsky, J. J. and White, T. J. (1990). PCR Protocols: A Guide to Methods and Applications. Academic Press, New York.

MSDS INFORMATION

Material Safety Data Sheets (MSDSs) are provided online at *http://www.genomics.agilent.com*. MSDS documents are not included with product shipments.